



DairyPal KNG

Rapid Test For The

Kanamycin, Neomycin, Gentamycin

In Milk



The Watar Bio DairyPal KNG KIT SC124 is a rapid test that detect simultaneously the presence of either Kanamycin Neomycin Gentamycin in a raw cow milk sample.



This DairyPal KNG KIT SC124 is specially designed to meet the European Union(EU) and Russian Federation(RF) regulation.

ADVANTAGES



MULTIPLE

Screening of 3 kinds of antibiotics in one single test



QUICK

9min per test



SENSITIVE

Detection limit more sensitive than MRLs



WIDELY USED

Adapt to EU and RF regulations



RELIABLE & STABLE

Performance robust and stable



USER FRIENDLY

Operate easily on site, in truck or lab




COST EFFECTIVELY

Save cost, time, labor

HOW TO OPERATE



Add milk

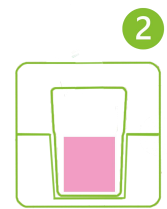

Incubate 3min at 45°C



Insert the Test



Incubate 6min at 45°C



Read Result



DETECTION LIMIT


ANTIBIOTICS	DETECTION LIMIT (PPB)
Kanamycin	10
Neomycin	50
Gentamycin	10

INSTRUMENT SUGGESTION

DRY INCUBATOR DairyPal Mini-B



READ THE TEST

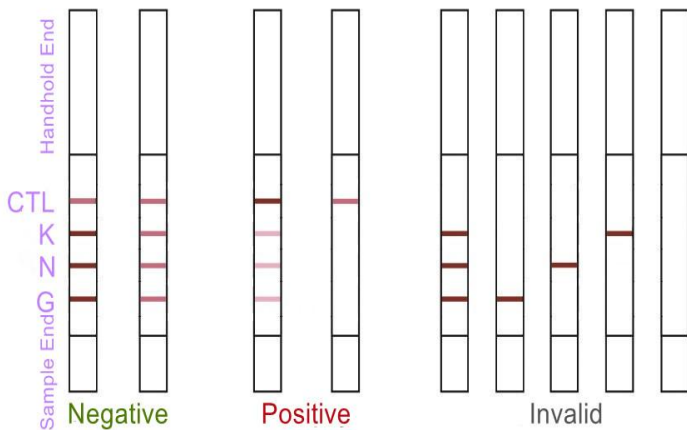

TEST READER DairyPal TSR-100



WatarBio

No. 12,
Hongzhuang south Road,
Yanta District, Xi'an,
China

+86-13679259477
info@watarbio.com
www.watarbio.com